Antibody Separation and Analysis





Better Surface Chemistry for Better Separation

Sepax Technologies, Inc.

Sepax Technologies, Inc. develops and manufactures products in the area of chemical and biological separations, bio-surfaces and proteomics. Sepax product portfolio includes 1) liquid chromatography columns and media, 2) SPE and Flash chromatography columns and tubes, 3) bulk resin for preparative separation and process chromatography, and 4) natural product and Chinese traditional medicine separation and purification.

A leader in Biological Separations

Sepax develops and manufactures wide range of biological separation products using both silica and polymeric resins as the support. The selection of particle size is from 1 μ m to 100 μ m and pore size from non-porous to 2000 Å. Unique and proprietary resin synthesis and surface technologies have been developed for solving the separation challenges in biological area.

Bioseparation Products

Size Exclusion SRT[®] SRT[®]-C Nanofilm[®] ZenixTM ZenixTM-C <u>Ion-exchange</u> Proteomix[®] Glycomix[™] <u>Antibody Separation</u> AntibodixTM Carbohydrate Separa

Carbohydrate Separation Carbomix[®]

Analytical, Semi-prep and Preparative







AntibodixTM NP Phases

General Description

Antibodix NP columns are specially designed for high resolution, high efficiency, and high recovery separations of antibodies. The packing support is composed of a rigid, spherical, highly cross-linked poly (styrene divinylbenzene) (PS/DVB) non-porous bead. The non-porous resin has particle size of 1.7, 3, 5 and 10 µm. The PS/DVB resin surface is grafted with a highly hydrophilic, neutral polymer thin layer with the thickness in the range of nanometer. On the top of the hydrophilic layer, weak cation-exchange functional groups are attached via a proprietary chemistry, resulting in high capacity ion-exchange layer.

Chemical Structure of Antibodix Resins

The chemical structure of Antibodix NP phases is composed of a rigid PS/DVB core, a densely packed, nanometer thick, hydrophilic coating, and a uniform weak cation exchange layer, as shown in Figure 1.







Highlights of Antibodix NP Resins

- · High separation efficiency and resolution
- Particle size selection of 1.7, 3, 5 and 10 μm
- Mono-dispersed particles
- Medium capacity
- High pressure tolerance: 4,000, 6,000, 8,000 and 12,000 psi for 10, 5, 3 and 1.7 µm resins, respectively
- Wide pH range: 2-12
- · High resolving power for slightly differed structures of monoclonal antibodies
- 1.7 and 3 μm particles are best suitable for high efficiency separation of proteins and MAbs

· Suitable for both analytical and scale-up separations of monoconal antibodies and other proteins

High Separation Efficiency

Antibodix NP resins have three unique features: non-porous particle, hydrophilic surface and a uniform layer of ionexchange functional groups, which enables high efficiency separations. Figure 2 is an example for separation of three proteins: ribonuclease A, cytochrome C, and lysozyme by Antibodix NP5 column. The average efficiency of three proteins reaches 132,000 of plates.



Figure 2. Separation of a protein mixture by Antibodix NP5 phase

Lot-to-Lot Reproducibility

With well-controlled resin production and the surface chemistry, manufacturing of Antibodix NP resins is highly reproducible. The typical variation of the retention time is less than 5% from batch to batch. One example is shown in Figure 3 for the production of three lots of Antibodix NP5 resins.

Chemical Dimension Availability

The column dimensions of Antibodix NP products are 0.75, 2.1, 3.0, 4.6, 7.8, 10, and 21.2 mm l.D., and 2, 3, 5, 10, 15, 25, and 30 cm length. We also offer custom-made columns.







Technical Specifications

Products	Particle size (µm)	Pressure limit (psi)	pH range	Temperature limit (⁰ C)	DBC* (mg/mL)
Antibodix-NP1.7	1.7	12,000	2-12	80	26.53 <u>+</u> 0.41
Antibodix-NP3	3	8,000	2-12	80	19.50 <u>+</u> 0.74
Antibodix-NP5	5	6,000	2-12	80	13.41 <u>+</u> 0.17
Antibodix-NP10	10	4,000	2-12	80	8.41 <u>+</u> 0.40

Dynamic binding capacity tests conditions: Sample: 3.0 mg/mL Lysozyme in 20 mM sodium phosphate buffer, pH 6.0; Flow rate: 0.5 mL/min (0.25 mL/min for 1.7 µm); Detection: UV 254 nm. Test on 5 different resin lots.

Applications

Separation & Analysis
Monoclonal antibodies (MAb)
MAb derivatives
Modified MAb molecules
Other proteins and peptides



Separation method development of a commercial monoclonal antibody sample

For a commercial monoclonal antibody sample, the separation conditions are critical for achieving optimized resolution. The key parameters of the separation conditions include salt concentration, pH and salt gradient. Figure 4 shows the separation of a commercial MAb sample, MAb-X22, with the mobile phase of 50 mM phosphate buffer, pH 6.0 at various gradients. Apparently the resolution is poor under those separation conditions.





0.0 1.0 2.0 3.0 4.0 5.0 6.0 7.0 8.0 9.0 10.0 11.0 12.0 13.0 14.0 15.0 16.0

Columns:	Antibodix-NP10 (10 μm, 4.6x250 mm)
Mobile phase:	A, 50 mM phosphate, pH 6.0
	B, A + 0.25 M NaCl
Gradient:	A) 0-100% B in 30 min
	B) 0-100%B in 45 min
	C) 0-30% B in 30 min
Flow rate:	0.8 mL/min
Sample:	MAb-X22
Injection:	10 μL (1.5 mg/mL)
Temperature:	25 °C
Detection:	UV 214 nm

After we optimized the separation conditions, the resolution of MAb-X22 is much better than that in Figure 4, as shown in Figure 5. Further on, we investigated the impact of gradient from shallow to deeper (25% to 60%B for 30 min). The trend is that more resolution comes with shallower gradient. However, the retention time increases when the gradient becomes shallower.

The initial salt concentration has great impact on the resolution of MAb samples. Figure 6 shows the separations of MAb-X22 sample with initial salt concentrations at 5, 10 and 20 mM phosphate at pH 7.5. 20 mM phosphate salt resulted in poor resolution, indicating that the initial salt concentration is very sensitive for resolving the fine structures of the MAb samples.

Figure 5. Separation of MAb-X22 with optimized conditions



Figure 6. The Impact of initial salt on the separation of MAb-X22



Columns:	Antibodix-NP10 (10 µm, 4.6x250 mm)				
Mobile phase:	A, Phosphate buffer, pH 7.5				
	B, A + 0.1 M NaCl				
Initial salt:	A/B/C=20/10/5 mM phosphate				
Gradient:	15-65%B in 60 min				
Flow rate:	0.8 mL/min				
Sample:	MAb-X22				
Injection:	10 μL (1.5 mg/mL)				
Temperature:	25 °C				
Detection:	UV 214 nm				

Figure 7 shows the impact of pH on the separation of MAbs. When pH decreased from 7.5 to 7.0, the resolution of the basic compound from the main component had some improvement with the compromise of longer retention time.





Figure 8. Separation of MAb-X22 with various gradients at pH 7



5

Figure 9 presents a separation profile of the monoclonal antibody. The main peak has a plate number of 12,170.





Quality control of a peptide drug molecule

Figure 10. Analysis of a commercial peptide drug molecule



Product Information for Antibodix NP Phase

ID x Length (mm)	P/N	Column Material	Phase	ID x Length (mm)	P/N	Column Material
7.8 x 75 7.8 x 50 4.6 x 100 4.6 x 50 4.0 x 10 (Guard) 2.1 x 50 2.0 x 10 (Guard) 4.6 x 50 Precolumn Filter	602NP2-7807 602NP2-7805 602NP2-4610 602NP2-4605 602NP2-4001C 602NP2-2105 602NP2-2001C 602NP2-2001C 602NP2P-4605 102000-P356	SS SS/PEEK SS/PEEK SS/PEEK SS/PEEK SS/PEEK SS/PEEK	Antibodix NP3 (3 μm)	7.8 x 75 7.8 x 50 4.6 x 150 4.0 x 10 (Guard) 2.1 x 50 2.0 x 10 (Guard) 4.6 x 50 Precolumn Filter	602NP3-7875 602NP3-7805 602NP3-4615 602NP3-4605 602NP3-4001C 602NP3-2105 602NP3-2001C 602NP3-2001C 602NP3P-4605 102000-P356	SS SS/PEEK SS/PEEK SS/PEEK SS/PEEK SS/PEEK SS/PEEK
10.0×250 7.8 x 150 7.8 x 75 4.6 x 250 4.6 x 50 4.0 x 10 (Guard) 2.1 x 150 2.1 x 50 2.0 x 10 (Guard) 4.6 x 250 4.6 x 50 Precolumn Filter	602NP5-10025 602NP5-7815 602NP5-7807 602NP5-4625 602NP5-4605 602NP5-4001C 602NP5-2115 602NP5-2105 602NP5-2001C 602NP5P-4625 602NP5P-4605 102000-P355 ep and preparative c 602NP5-21225	SS SS SS/PEEK SS/PEEK SS/PEEK SS/PEEK SS/PEEK SS/PEEK SS/PEEK SS/PEEK	Antibodix NP10 (10 μm)	10.0×250 7.8 x 150 7.8 x 75 4.6 x 250 4.6 x 50 4.0 x 10 (Guard) 2.1 x 250 2.1 x 50 2.0 x 10 (Guard) 4.6 x 250 4.6 x 50 Precolumn Filter Semi- 21.2 x 250 24.0 x 450	602NP10-10025 602NP10-7815 602NP10-7807 602NP10-4625 602NP10-4605 602NP10-4001C 602NP10-2115 602NP10-2105 602NP10-2001C 602NP10P-4625 602NP10P-4605 102000-P355 prep and preparative co 602NP10-21225	SS SS SS/PEEK SS/PEEK SS/PEEK SS/PEEK SS/PEEK SS/PEEK SS/PEEK SS/PEEK
- - - - - - - - - - - - - - - - - - -	7.8 x 150 7.8 x 75 4.6 x 250 4.6 x 50 4.0 x 10 (Guard) 2.1 x 150 2.1 x 150 2.1 x 50 2.0 x 10 (Guard) 4.6 x 250 4.6 x 50 Precolumn Filter 21.2 x 250 21.2 x 150	7.8 x 150 602NP5-7815 7.8 x 75 602NP5-7807 4.6 x 250 602NP5-4625 4.6 x 50 602NP5-4605 4.0 x 10 602NP5-4001C (Guard) 602NP5-2115 2.1 x 150 602NP5-2105 2.0 x 10 602NP5-2001C (Guard) 602NP5-2001C 4.6 x 250 602NP5P-4625 4.6 x 50 602NP5P-4605 Yrecolumn 102000-P355 Filter Semi-prep and preparative c 21.2 x 250 602NP5-2125	7.8 x 150 602NP5-7815 SS 7.8 x 75 602NP5-7807 SS 4.6 x 250 602NP5-4625 SS/PEEK 4.6 x 50 602NP5-4605 SS/PEEK 4.6 x 50 602NP5-4001C SS/PEEK 4.0 x 10 602NP5-4001C SS/PEEK 2.1 x 150 602NP5-2115 SS/PEEK 2.1 x 150 602NP5-2105 SS/PEEK 2.1 x 50 602NP5-2001C SS/PEEK 2.0 x 10 602NP5-2001C SS/PEEK (Guard) 602NP5P-4625 SS/PEEK 4.6 x 250 602NP5P-4625 SS/PEEK 4.6 x 50 602NP5P-4605 SS/PEEK Yrecolumn Filter 102000-P355 SS/PEEK 21.2 x 250 602NP5-21225 SS 21.2 x 150 602NP5-21215 SS	7.8 x 150 602NP5-7815 SS 7.8 x 75 602NP5-7807 SS 4.6 x 250 602NP5-4625 SS/PEEK 4.6 x 50 602NP5-4605 SS/PEEK 4.6 x 50 602NP5-4001C SS/PEEK 4.0 x 10 602NP5-4001C SS/PEEK (Guard) 602NP5-2115 SS/PEEK 2.1 x 150 602NP5-2105 SS/PEEK 2.1 x 50 602NP5-2105 SS/PEEK 2.1 x 50 602NP5-2001C SS/PEEK 2.0 x 10 602NP5-2001C SS/PEEK (Guard) 602NP5-4625 SS/PEEK 4.6 x 250 602NP5P-4625 SS/PEEK Yrecolumn 102000-P355 SS/PEEK Filter 102000-P355 SS/PEEK Semi-prep and preparative columns SS 21.2 x 250 602NP5-21225 SS 21.2 x 150 602NP5-21215 SS	7.8 x 150 602NP5-7815 SS 7.8 x 150 7.8 x 75 602NP5-7807 SS 7.8 x 75 4.6 x 250 602NP5-4625 SS/PEEK 4.6 x 250 4.6 x 50 602NP5-4605 SS/PEEK 4.6 x 50 4.0 x 10 602NP5-4001C SS/PEEK 4.6 x 50 2.1 x 150 602NP5-2115 SS/PEEK 4.0 x 10 2.1 x 50 602NP5-2105 SS/PEEK 2.1 x 250 2.1 x 50 602NP5-2001C SS/PEEK 2.1 x 50 2.0 x 10 602NP5-2001C SS/PEEK NP10 2.0 x 10 (Guard) 602NP5P-4625 SS/PEEK 4.6 x 250 4.6 x 250 4.6 x 50 602NP5P-4625 SS/PEEK 4.6 x 50 Yerecolumn Filter 102000-P355 SS/PEEK 4.6 x 50 Precolumn Filter 102000-P355 SS/PEEK 21.2 x 250 21.2 x 250 21.2 x 250 602NP5-21225 SS 21.2 x 250 21.2 x 150	7.8 x 150 602NP5-7815 SS 7.8 x 75 602NP5-7807 SS 4.6 x 250 602NP5-4625 SS/PEEK 4.6 x 250 602NP5-4605 SS/PEEK 4.6 x 50 602NP5-4605 SS/PEEK 4.6 x 50 602NP5-4001C SS/PEEK 4.6 x 50 602NP5-2115 SS/PEEK 2.1 x 150 602NP5-2105 SS/PEEK 2.1 x 150 602NP5-2001C SS/PEEK 2.0 x 10 602NP5-2001C SS/PEEK (Guard) 602NP5P-4625 SS/PEEK 2.0 x 10 602NP5-2001C SS/PEEK (Guard) 602NP5P-4625 SS/PEEK 4.6 x 250 602NP5P-4625 SS/PEEK 4.6 x 50 602NP5P-4625 SS/PEEK 4.6 x 50 602NP5P-4605 SS/PEEK 4.6 x 50 602NP5P-4605 SS/PEEK Precolumn 102000-P355 SS/PEEK Precolumn Filter 102000-P355 S2/PEEK SS 21.2 x 250 602NP10-2125 21.2 x 250 602NP5-2125 SS 21.2 x 250 602NP10-21225

* SS means Stainless steel.

** Precolumn Filters comes with 0.5 μm PEEK frit for 102000-P356 and 2.0 μm PEEK frit for 102000-P355.

*** Other column dimensions and custom-made column dimensions are available.



Precolumn Filter





How to Order

It's fast and easy to order from the Sepax on-line store at: www.sepax-tech.com

Or contact Sepax Sales Department by **Phone:** (302) 366-1101 1-877-SEPAX-US **Fax:** (302) 366-1151

E-mail: sales@sepax-tech.com

Mail:

5 Innovation Way, Delaware Technology Park Newark, Delaware 19711 USA

Sepax Technologies, Inc. 5 Innovation Way, Newark, Delaware 19711, USA Tel: (302) 366-1101 Fax: (302) 366-1151 www.sepax-tech.com